




Icon-NMR Manual

(updated 2025.02.21)



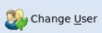

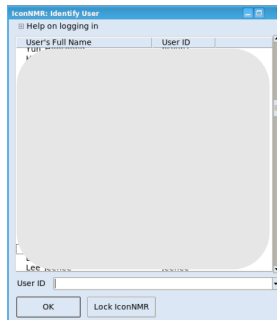

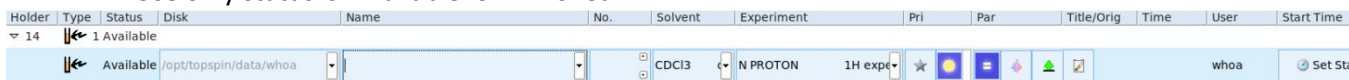

- Put NMR sample on tray
 - Remember holder number
- Login
 - Click  if login window is not appeared





- Double click available holder number
 - Use only status is 'Available' or 'Finished'



- Fill out below information

- Name : dataset name // Alphabet, number, hyphen, underbar only
- No. : experiment number
- Solvent : choose solvent
- Experiment : read description next to experiment name
- Priorty :  ignore star icon
- Day/Night :  Day (09:00 ~ 24:00).  Night (00:00 ~ 09:00)
 - Check rule at nmrchem.kaist.ac.kr [Policy]
- Parameter : modify parameter if you know what is
 - NS : Number of scan, D1 : Relaxation delay
 - AQ, O1P, 1SW
 - Before modification , after modification 
- Lock/Shim/Tune: modify if you need 
- Print : ignore 
- Title : Comment 

Status	설명	Can delete by other user
Available	사용가능	O
Queued	예약됨	X
Running	분석중	X
Finished	분석완료	O
Failed	분석실패	X

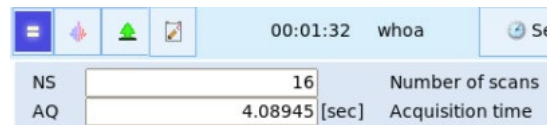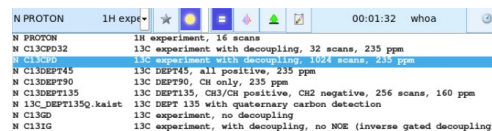


- Click 'Submit' icon

- Time : expected time
- Start time : expected starting time
- Submit : submit queue . Cancel :cancel queue, Edit : edit parameters
- Add : add additional experiments at one holder number
- Copy : make new queues with same parameter setting



00:00:51 whoa 21:29 Wed Apr 06 2022

- Click 'Change User'

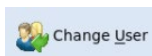

Report quickly if LED is red